



**BMA BIOMEDICALS**

BMA BIOMEDICALS  
Rheinstrasse 28-32  
CH-4302 Augst (Switzerland)  
Phone: ++41 61 811 6222  
Fax: ++41 61 811 6006  
info@bma.ch  
www.bma.ch

---

## **Staining Procedure For Frozen Tissue Sections With Product T-1056**

### **Protocol with frozen, ice-cold acetone-fixed sections:**

The whole procedure is performed at room temperature

1. Wash in PBS
2. Block endogenous peroxidase
3. Wash in PBS
4. Block with 10% normal goat serum in PBS for 30min. in a humid chamber
5. Incubate with primary antibody (dilution see datasheet) for 1h in a humid chamber
6. Wash in PBS
7. Incubate with secondary antibody (peroxidase-conjugated goat anti mouse IgG+IgM (H+L) minimal-cross reaction to human) for 1h in a humid chamber
8. Wash in PBS
9. Incubate with AEC substrate (3-amino-9-ethylcarbazol) for 12min.
10. Wash in PBS
11. Counterstain with Mayer's hemalum

For further information and details see *technical information*