



BMA BIOMEDICALS

BMA BIOMEDICALS
Rheinstrasse 28-32
CH-4302 Augst (Switzerland)
Phone: ++41 61 811 6222
Fax: ++41 61 811 6006
info@bma.ch
www.bma.ch

Staining Procedure For Frozen And Paraffin-Embedded Tissue Sections With Product T-1016

Protocol with frozen, ice-cold acetone-fixed sections:

The whole procedure is performed at room temperature

1. Wash in PBS
2. Block endogenous peroxidase
3. Wash in PBS
4. Block with 10% normal goat serum in PBS for 30min. in a humid chamber
5. Incubate with primary antibody (dilution see datasheet) for 1h in a humid chamber
6. Wash in PBS
7. Incubate with secondary antibody (peroxidase-conjugated goat anti mouse IgG+IgM (H+L) minimal-cross reaction to human) for 1h in a humid chamber
8. Wash in PBS
9. Incubate with AEC substrate (3-amino-9-ethylcarbazol) for 12min.
10. Wash in PBS
11. Counterstain with Mayer's hemalum

Protocol with formalin-fixed, paraffin-embedded sections:

The whole procedure is performed at room temperature

1. Deparaffinize and rehydrate tissue section
2. Incubate the tissue section with proteinase K for 7min.
3. Wash in distilled water
4. Block endogenous peroxidase
5. Wash in PBS
6. Block with 10% normal goat serum in PBS for 30min. in a humid chamber
7. Incubate with primary antibody (dilution see datasheet) for 1h in a humid chamber
8. Wash in PBS
9. Incubate with secondary antibody (peroxidase-conjugated goat anti mouse IgG+IgM (H+L) minimal-cross reaction to human) for 1h in a humid chamber
10. Wash in PBS
11. Incubate with AEC substrate (3-amino-9-ethylcarbazol) for 12min.
12. Wash in PBS
13. Counterstain with Mayer's hemalum

For further information and details see ***technical information***