



BMA BIOMEDICALS

BMA BIOMEDICALS
Rheinstrasse 28-32
CH-4302 Augst (Switzerland)
Phone: ++41 61 811 6222
Fax: ++41 61 811 6006
info@bma.ch
www.bma.ch

Staining Procedure For Frozen Mouse Tissue Sections With Product T-2109

Protocol with frozen, ice-cold acetone-fixed sections:

The whole procedure is performed at room temperature

1. Wash in PBS
2. Block endogenous peroxidase
3. Wash in PBS
4. Block with 10% normal goat serum in PBS for 30min. in a humid chamber
5. Incubate with primary antibody (dilution see datasheet) for 1h in a humid chamber
6. Wash in PBS
7. Incubate with secondary antibody (peroxidase-conjugated goat anti rat IgG (H+L) minimal-cross reaction to mouse) for 1h in a humid chamber
8. Wash in PBS
9. Incubate with AEC substrate (3-amino-9-ethylcarbazol) for 12min.
10. Wash in PBS
11. Counterstain with Mayer's hemalum

For further information and details see *technical information*