



BMA BIOMEDICALS



Peninsula Laboratories

---

## Goat anti Rabbit IgG Antiserum, pretitrated

This antiserum has been titrated against normal rabbit serum and is designed to be used specifically in conjunction with pretitrated Normal Rabbit Serum (catalog #T-4577) for precipitation of "bound" counts in the titration of antisera or in RIAs.

---

### TECHNICAL AND ANALYTICAL CHARACTERISTICS

**Product number:** T-4576 (Lot A19PO2158)

**Host species:** Rabbit IgG

**Format:** Serum, lyophilized, packaged under nitrogen

#### **Reconstitution Information:**

Rehydrate the lyophilized powder with 50 ml of assay buffer.

#### **Recommended Use:**

- Add 100µl of rehydrated Goat anti-Rabbit IgG to each tube
- Add to each tube, 100µl of pretitrated Normal Rabbit Serum (catalog #T-4577), and vortex
- Allow the precipitate to form for 2 hours at room temperature (or 6 hours at 4°C)
- Centrifuge the tube at 1700xg for 20 min to collect the precipitate
- Collect the precipitate and carefully aspirate the supernatant
- Count the pellet in a scintillation well gamma counter

#### **Related Products:**

T-4577: Normal Rabbit Serum (pretitrated)  
Y-1000: SEP-Columns (200mg, 50 per box)  
Y-1040: Buffer A for sample extraction  
Y-1045: Buffer B for sample extraction  
Y-1050: RIA Buffer

---

This product contains no preservative and is intended for laboratory use and research purposes only. Not for diagnostic or therapeutic applications.

T-4576

07.08.2024