



BMA BIOMEDICALS



Peninsula Laboratories

T-5052

Normal Guinea Pig Serum (Pretitered)

This serum has been designed to be used specifically in conjunction with pretitered Goat anti-Guinea Pig IgG (GAGGG) for precipitation of "bound" counts in the titration of antisera or in RIA's.

Reconstitution Information:

Rehydrate the lyophilized powder with 50ml of buffer, this should be the same buffer used in the assays.

Recommended Use:

After the addition of 100µl of GAGGG to each tube, add 100µl of the rehydrated NGS to each tube and vortex the contents. Allow the precipitates to form for 2 hours at room temperature. Centrifuge the tubes at 1700xg for 20 minutes and then carefully aspirate off the supernatant. Count the pellet in a scintillation well gamma counter.

T-5052

17.3.2021